

MEMBRAN FILTRATION KIT

M- NUTRIENT TTC BROTH 3,6 CYROVIAL

INTENDED USE:

Nutrient TTC Broth is used for the general cultivation of less fastidious microorganisms. Membrane filter technique is more convenient to work with high volume samples and count as well as conventional procedures.

PRINCIPLE AND INTERPRETATION:

Membrane filter technique is an effective, accepted technique for testing fluid samples for microbiological contamination. It involves less preparation than many traditional methods, and is one of a few methods that will allow the isolation and enumeration of microorganisms.

TEST PROCEDURE:

Sample Volume : A sample volume of 50 to 100 ml should be selected.

Filtration Technique :

1. The filter set body is sterilized in an autoclave at 121 oC for 15-30 minutes and the filter assembly is installed.
2. Membrane filter is taken from the sterile pack with a sterile clamp.
3. Carefully placed in the container with the checkered side of the filter on top.
4. Turn on the vacuum and allow the sample to draw completely through the filter.
5. While the filtration is in progress, the 60 mm petri dish with sterile pad opens.
6. 3.6 ml of m-Endo Broth are poured into the petri dish.
7. After the filtration process is finished, the membrane filter is taken carefully with the help of a sterile pliers.
8. Place the membrane filter into the prepared Petri dish.
9. Incubate at the proper temperature and for the appropriate time period.

COMPOSITION OF BROTH:

Ingredients	Gr/Liter
Peptone	5 gr
Sodium chloride	5 gr
Peptone-B	1,5 gr
Yeast extract	1,5 gr

***Formula adjusted, standardized to suit performance parameters

pH: 7,4 ± 0,2

QUALITY CONTROL OF MEDIA:

1.Sterility Control:

Incubation 48 hours at 30-35°C and 72 hours at 20-25°C: NO GROWTH

2.Physical/Chemical Control

pH: 7,4 ± 0,2

Apperance: Light amber

3.Microbiological Control: Incubation at 35± 2 °C during 24 h

Microorganism	Inoculum (CFU)	Results	
		Growth	Reaction
Pseudomonas aeruginosa ATCC 27853	10-100	Good	Reddish colonies
Enterococcus faecalis ATCC 29212	10-100	Good	Reddish colonies
E.coli ATCC 25922	10-100	Good	Reddish colonies
Salmonella typhimurium ATCC 14028	10-100	Good	Reddish colonies

PRECAUTIONS:

RTA.KK.670 Revision Date/Revision Number:-/0 Issue Date: 15.03.2016

For professional use only. Do not use plates if they show evidence of microbial contamination, discoloration, drying, cracking or other signs of deterioration.

STORAGE CONDITIONS AND SHELF LIFE:

Store the prepared medium at 2 - 12°C. Use before expiry date on the label. Do not use beyond stated expiry date.

DISPOSAL:

Incubated prepared medium may contain active bacteria and micro-organisms. Do not open infected medium. Infected plate should be autoclaved, incinerated or opened and soaked in a chlorine-based disinfectant (liquid bleach) for 20 minutes prior to disposal.

PACKAGING:

Katalog Number: 06253

Packaging:Box

Content: 100 cyrovial, 100 membran filters and petri dishes with pad/each box

REFERENCES:

1. Fifiield, C. W., and C. P. Schaufus. 1958. J. Am. Water Works Assoc. 50:193-196.
2. Eaton, A. D., L. S. Clesceri, and A. E. Greenberg (eds.). 1998. Standard methods for the examination of water and wastewater, 20th ed. American Public Health Association, Washington, D.C.
3. Downes, F. P., and K. Ito (eds.). 2001. Compendium of methods for the microbiological examination of foods, 4th ed. American Public Health Association, Washington, D.C.
4. Bordner, R., and J. Winter (eds.). 1978. Microbiological methods for monitoring the environment, water, wastes. EPA-600/8-78-017. Environmental Monitoring and Support Laboratory, Office of Research and Development, U. S. Environmental Protection Agency, Cincinnati, OH.
5. U. S. Environmental Protection Agency. 1992. Manual for the certification of laboratories analyzing drinking water. EPA-814B92-002. Office of Ground Water and Technical Support Division, U. S. Environmental Protection Agency, Cincinnati, OH.



Aseptic Sterile



Batch Code



Catalogue Number



Negative Controls



Positive Controls



Use by



Temperature Limitation



Do not reuse



Contains sufficient for <n> tests



Look at user manual



Manufacturer